

CI-hAELVi

Cat. No.: INS-CI-1015



Biosafety Level
Level 1



Storage
vapor phase of liquid nitrogen



Culture Conditions
huAEC Coating Solution (INS-SU-1018)
huAEC Medium (INS-ME-1013)

General Information

CI-hAELVi is a human alveolar epithelial cell line that was immortalized using the [CI-SCREEN](#) technology ([Lipps et al. 2018; Nat Comm](#)) and described by [Kuehn et al. 2016; ALTEX](#). The cell line exhibits Type-I-like properties (expression of caveolin-1, absence of surfactant protein C) and expresses epithelial markers (occluding, ZO-1). For a list of selected references refer to page 3.

Organism: Homo sapiens (human)

Tissue: lung

Growth properties: adherent

Cell culture media and reagents

Product	Cat. No.	Volume
huAEC Medium (includes basal medium and supplements)	INS-ME-1013	500ml
huAEC Coating Solution	INS-SU-1018-20ml INS-SU-1018-100ml	20ml 2x50ml
Freezing medium	INS-SU-1004	30ml

Note: The medium does not contain antibiotics. However, it may simply be supplemented with standard antibiotics.

Intended Use

This product is intended for in vitro research use only. It is not intended for any animal or human therapeutic or diagnostic use.

Quality control

Each vial contains $\geq 5 \times 10^5$ cells. Viability is $\geq 80\%$. Cells are negative for mycoplasma contamination. The source material is tested negative for HIV, HBV and HCV.

Upon arrival

Cells are routinely shipped on dry ice. Check all containers for leakage and breakage. Check if cells arrived frozen. After arrival, store the cryopreserved cells in liquid nitrogen vapor, or seed them immediately (please see page 2).

Note: Cells may be stored at -80°C for short periods (< 2 days), but this results in reduced viability and irreversible cell damage.

Coating of cell culture vessels

Material:

- Cell culture vessel(s)
- PBS
- huAEC Coating Solution (INS-SU-1018)

Protocol:

- 1) Cover the cell culture vessel with coating solution (see table below for the required volume).
- 2) Incubate the cell culture vessel for at least 2h (up to overnight) at 37°C in the incubator.
- 3) Aspirate coating solution.
- 4) Wash once with PBS and aspirate PBS.

Note: Coated cell culture vessels may be stored sealed at $2-8^\circ\text{C}$ for up to 7 days.

Vessel	Surface area (cm ²)	Volume (ml)
T75	75	2.5
T25	25	1.4
6well	9.6	0.7
12well	3.5	0.25
24 well	1.9	0.1
96well	0.32	0.05

Medium storage and preparation

Medium including supplements is shipped cooled at 4-8°C. Store reagents according to the instructions below upon arrival.

Material:

- huAEC Medium INS-ME-1013. Includes Basal Medium (500ml) and Supplements (30ml; INS-ME-1013BS).

Storage:

- Store Basal Medium at 4-8°C
- Store Supplements at -20°C
- Store completed medium (Basal Medium plus Supplements) at 4-8°C. Completed medium is stable for at least 1 month at 4-8°C.

Protocol:

- 1) Thaw Supplements at 15-25°C.
- 2) Add 30ml Supplements to 500ml Basal Medium and store at 4-8°C. Completed medium is stable for at least 1 month at 4-8°C.

Note: Supplements may be aliquoted and stored at -20°C before completing the medium. For example, aliquot 5×6ml and then add 6ml Supplements to 100ml Basal Medium.

Recover cryopreserved cells

Do not thaw the cells until the recommended medium and coated flasks are on hand. For initial recovery (after delivery), we recommend thawing the cells on a T25 flask.

Material:

- huAEC solution (INS-SU-1018) coated cell culture vessels
- pre-warmed (37°C) complete medium (Basal Medium plus Supplements)
- 15ml tube

Protocol:

- 1) Add 4ml pre-warmed medium to a 15ml tube.
- 2) Quickly thaw the cryovial at 37°C in a water bath until only a few ice crystals are visible. Disinfect vial briefly with 70% Ethanol.
- 3) Transfer thawed cell suspension to the 15ml tube containing 4ml medium. Avoid excessive pipetting up and down.
- 4) Centrifuge cells at 200×g for 4min.
- 5) Aspirate supernatant.
- 6) Gently resuspend the cell pellet in complete medium.
- 7) Transfer cells in coated cell culture vessel and place in the incubator (37°C, 5% CO₂).
- 8) Change the medium after 2 days.

Routine Subculture

Change medium every 2 days and split the cells at 70-90% confluence. The split ratio after recovery from cryopreservation should not exceed 1:2. For routine maintenance, split ratio can be increased to 1:3 to 1:4.

Material:

- huAEC solution (INS-SU-1018) coated cell culture vessels
- pre-warmed (37°C) complete medium (Basal Medium plus Supplements)
- PBS
- Trypsin/EDTA solution (TE)

Protocol:

- 1) Aspirate medium.
- 2) Wash with PBS and aspirate PBS.
- 3) Add Trypsin/EDTA (TE) solution to the cells and incubate at room temperature or 37°C for 5-10min, or until the cells detach.
- 4) Examine the cells under a microscope. When the cells start to detach, gently tap the side of the vessel to loosen the remaining cells.
- 5) Resuspend cells in complete medium thereby inactivating the Trypsin/EDTA (TE) solution.
- 6) Transfer an aliquot of the cell suspension to a new coated cell culture vessel containing fresh complete medium.
- 7) Incubate at 37°C and 5% CO₂.

Vessel	Medium or PBS (ml)	TE (ml)
T75 flask	8-10	2
T25 flask	4-5	1
6well	1.5-2	0.5
12well	1	0.2
24 well	0.5	0.1
96well	0.1	0.05

Cryopreservation

Cell should be grown to 90% confluence before cryopreservation. Avoid full confluence before cryopreservation.

Material:

- Freezing medium (INS-SU-1004)
- PBS
- Trypsin/EDTA solution (TE)
- 2% FBS in PBS
- 15ml tube
- cryovial(s)
- freezing container ("Mr. Frosty" or similar)

Protocol:

- 1) Aspirate medium.
- 2) Wash with PBS and aspirate PBS.
- 3) Add Trypsin/EDTA (TE) solution to the cells and incubate at room temperature or 37°C for 5-10min, or until the cells attach.
- 4) Examine the cells under a microscope. When the cells start to detach, gently tap the side of the vessel to loosen the remaining cells.
- 5) Resuspend cells in 2% FBS in PBS and transfer to a 15ml tube.
- 6) Centrifuge cells at 200×g for 4min.
- 7) Aspirate supernatant and gently resuspend cell pellet in Freezing medium (approx. 1×10^6 cells/ml).
- 8) Transfer cell suspension into cryovial(s) and place them into a freezing container ("Mr. Frosty" or similar).
- 9) Place the freezing container at -70 to -80°C for 16-24h.
- 10) Transfer cryovials to liquid nitrogen vapor for long-term-storage.

Cell Culture on Permeable Inserts

Cells can be cultured on cell culture inserts at the Liquid-Liquid interface (LLI) or Air-Liquid interface (ALI).

Material:

- huAEC Coating Solution (INS-SU-1018) coated cell culture inserts (e.g., SABEU cellQART®, Corning Transwell® or similar products)
- standard multiwell plate (to hold the cell culture inserts)
- pre-warmed (37°C) complete medium (Basal Medium plus Supplements)
- PBS

Protocol:

- 1) Add the appropriate amount of coating solution to the required number of inserts.

Insert	Coating Volume (mL)	Wash Volume (mL)
6 well	0.6	1
12 well	0.25	0.4
24 well	0.05	0.1
96 well	0.025	0.05

- 2) Incubate the cell culture inserts for at least 2h (up to overnight) at 37°C in the incubator.
- 3) Aspirate coating solution.
- 4) Rinse once with PBS and aspirate PBS.
Note: Coated cell culture inserts may be stored sealed at 2-8°C for up to 7 days.
- 5) Harvest and count cells according to routine subculture protocol.
- 6) Seed 3×10^5 cells/cm² onto pre-coated cell culture inserts, e.g., 1×10^5 cells onto one 24-well insert.
- 7) Change medium every 2 days. When changing the media, always aspirate from the basolateral side first, then the apical side. When replacing the media, add in the opposite order (apical then basolateral).
- 8) For ALI culture: air-lift cells after 2-3 days, but earliest when cells are confluent.

Contact Information

InSCREENeX GmbH
Inhoffenstr. 7
38124 Braunschweig
Germany

Email: info@inscreenex.com

Website: inscreenex.com

Phone: +49 531 6181 5080

Fax: +49 531 6181 5002

Selected References

hAELVi cell line generation and characterisation:

Kuehn, Anna et al. "Human alveolar epithelial cells expressing tight junctions to model the air-blood barrier." *ALTEX* vol. 33,3 (2016): 251-60. doi:10.14573/altex.1511131 [\[Link\]](#)

hAELVi co-culture system with THP-1 macrophages:

Kletting, Stephanie et al. "Co-culture of human alveolar epithelial (hAELVi) and macrophage (THP-1) cell lines." *ALTEX* vol. 35,2 (2018): 211-222. [\[Link\]](#)

hAELVi culture in microfluidic-based acini-on-chips:

Artzy-Schnirman, Arbel, et al. "Capturing the Onset of Bacterial Pulmonary Infection in Acini-On-Chips." *Advanced Biosystems* (2019). [\[Link\]](#)

In vitro exposure of hAELVi cells to halogenated gases:

Blondonnet, Raïko, et al. "In Vitro Method to Control Concentrations of Halogenated Gases in Cultured Alveolar Epithelial Cells." *JoVE* 140 (2018): e58554. [\[Link\]](#)

Nanoparticle toxicity assessment in vitro using hAELVi cells:

Leibrock, Lars, et al. "Nanoparticle induced barrier function assessment at liquid-liquid and air-liquid interface in novel human lung epithelia cell lines." *Toxicology Research* (2019). [\[Link\]](#)

Release characteristics of siRNA from nanocarriers:

Thanki, Kaushik et al. "Mechanistic profiling of the release kinetics of siRNA from lipidoid-polymer hybrid nanoparticles in vitro and in vivo after pulmonary administration." *Journal of controlled release*. vol. 310 (2019): 82-93. [\[Link\]](#)

Biocompatibility testing of carrier systems for pulmonary delivery of anti-infective:

Ho, Duy-Khiet, et al. "Polysaccharide Submicrocarrier for Improved Pulmonary Delivery of Poorly Soluble Anti-infective Ciprofloxacin: Preparation, Characterization, and Influence of Size on Cellular Uptake." *Mol pharmaceutics* 15.3 (2018): 1081-1096. [\[Link\]](#)